SEPAX TECH.

YOUR SEPARATION PARTNER

Sepax Technologies

Represented by

Analab

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

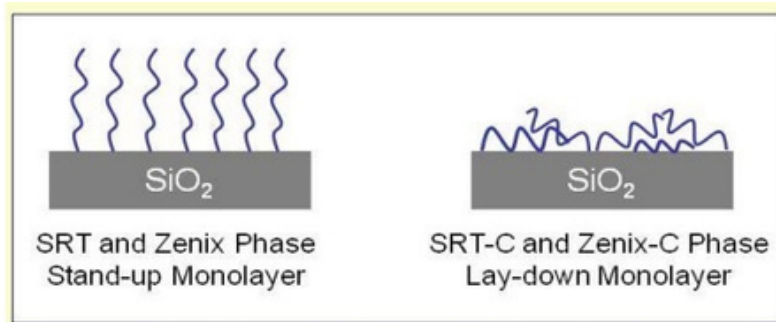
Characteristics	SRT	Zenix	SRT-C	Zenix-C
Particle size	5 μm	3 μm	5 μm	3 μm
Pore size (Å)	100, 150, 300, 500, 1000 & 2000	80, 100, 150, 300	100, 150, 300, 500, 1000 & 2000	80, 100, 150, 300
Resolution	High	Higher, short column for faster separation	High	Higher, short column for faster separation
Surface structure	Chemically bonded stand-up monolayer		Chemically bonded lay-down monolayer	
Recommended Sample Types	Monoclonal antibodies, proteins, peptides, nucleic acids, oligonucleotides, virus, and water-soluble polymers		"Tough samples" such as hydrophobic samples like insulin, membrane protein, antibody drug conjugates, proteins conjugated with hydrophobic molecules.	

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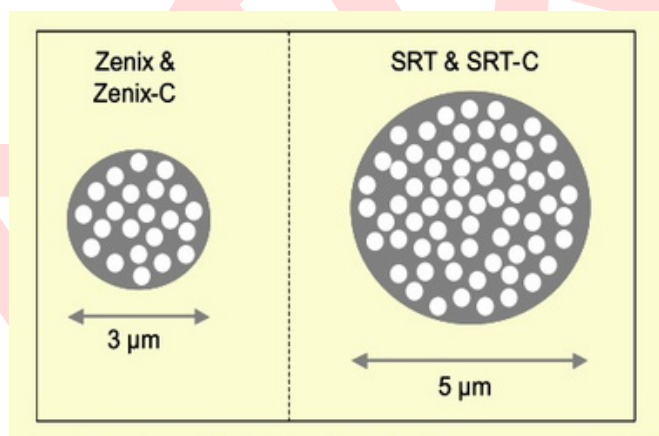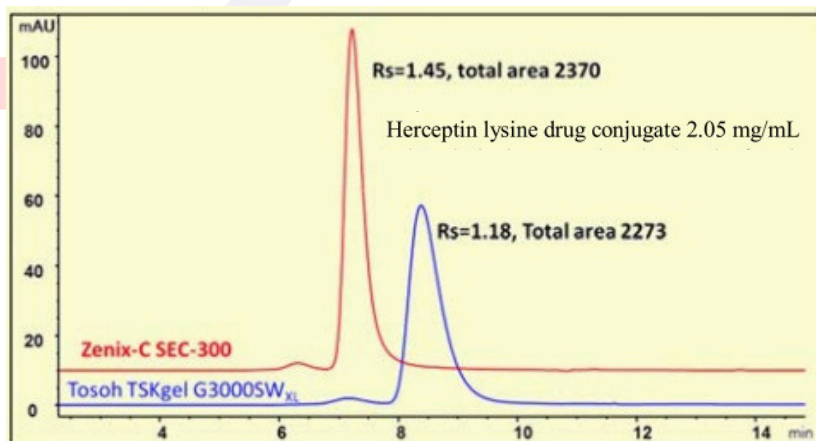Column: **Zenix - C SEC-300** (3 μm , 300 Å, 7.8 x 300 mm) & **Tosoh TSKgel G3000SWXL** (5 μm , 250 Å, 7.8 x 300 mm)
Mobile phase: 150 mM sodium phosphate buffer, pH 7.0 1 mL/min
Detection: UV 280 nm
Temperature: Ambient
Injection: 10 μL
Samples: Herceptin lysine drug conjugate 2.05 mg/mL
Disclaimer: TSKgel and Tosoh Bioscience are registered trademarks of Tosoh Corporation; Comparative separations may not be representative of all applications.

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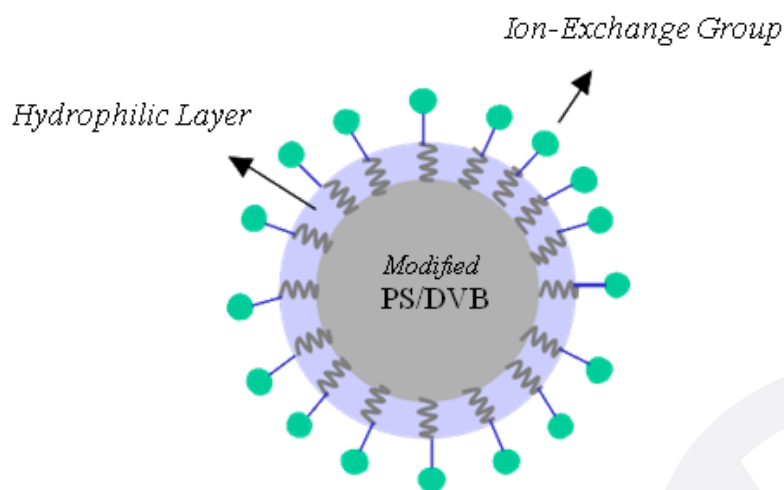
Phase	Proteomix [®] SCX, WCX
Dimensions	4.6 x 100 mm
Material	Non-porous PS/DVB beads grafted with a highly hydrophilic, neutral polymer thin layer.
Maximum temperature (°C)	~ 80
Mobile phase compatibility	Aqueous or a mixture of water and acetonitrile, acetone, or methanol

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