YOUR BIO SEPARATION PARTNER

Sepax Technologies

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Contact:

TEL : 02-2776-6931

MAIL : sales@analab.com.tw

LINE : @analab

SRT SEC-750

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
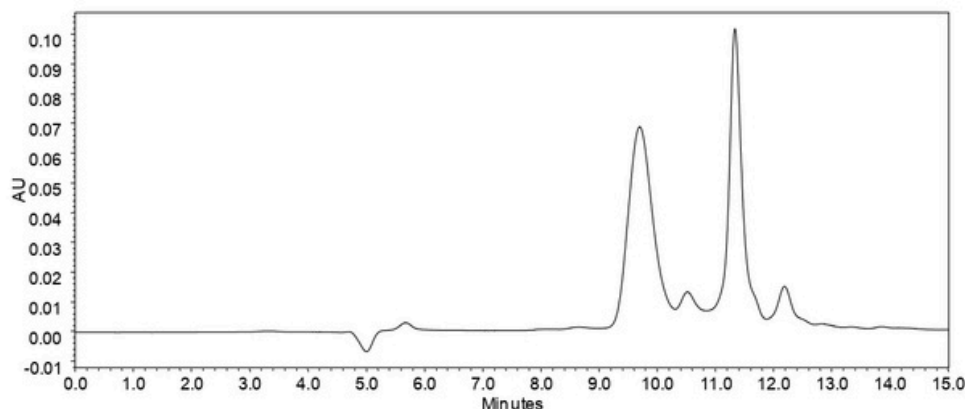
Separation of Thyroglobulin on SRT®-1000

Column: SRT®SEC-1000 (5 μm , 1000 \AA , 4.6 x 300 mm)
Mobile Phase: 150 mM phosphate buffer, pH 7
Flow Rate: 0.35 mL/min
Temperature: Ambient
Sample: rhIL-10 sR
Concentration: 1 mg/mL
Injection Volume: 3 μL

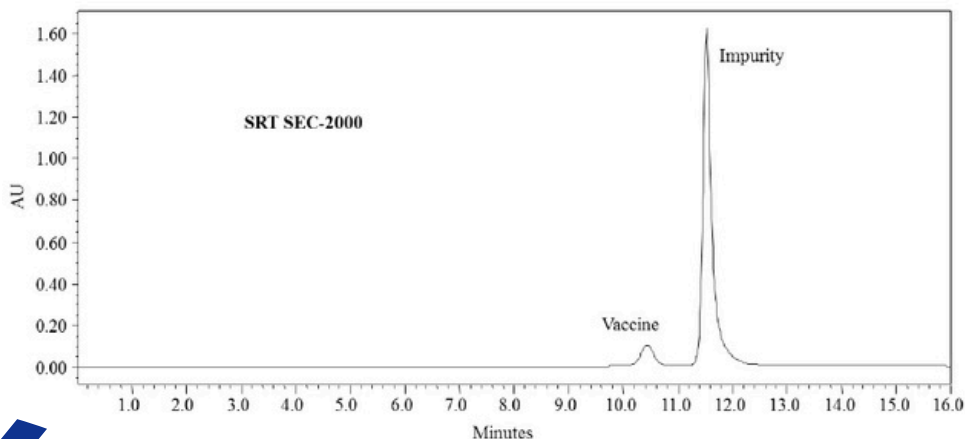Analysis of Vaccine Sample on SRT®SEC-2000

Column: SRT® SEC-2000
Mobile phase: 150 mM phosphate buffer, pH 7.0
Flow rate: 1.0 mL/min
Temperature: 25°C
Sample injection: 20ul
Sample: Vaccine
MW~80,000 KD

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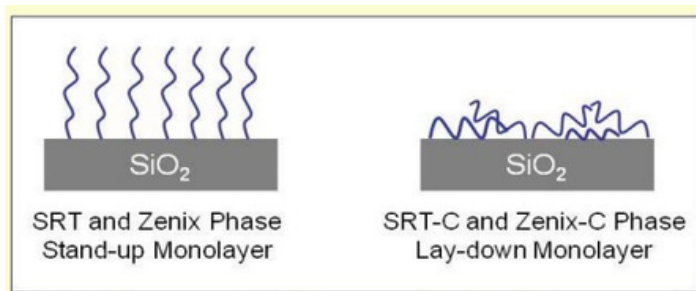
Ex: **hydrophobic proteins** like insulin, membrane protein, monoclonal antibodies derivatized with polymer branches, e.g. polypeptide, PEG.

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
Herceptin-Lysine Drug Conjugate Analysis

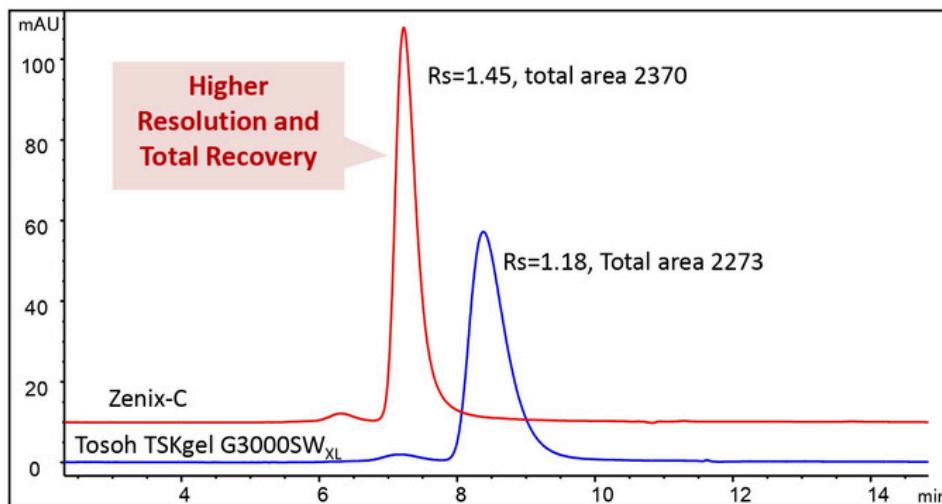Zenix-C SEC-300 vs. Tosoh TSKgel G3000SWXL

Column: Zenix SEC-300 (3 μ m, 300 \AA , 7.8 x 300 mm)

Mobile phase: 150 mM phosphate buffer, pH 7.0;

Flow rate: 1 mL/min;

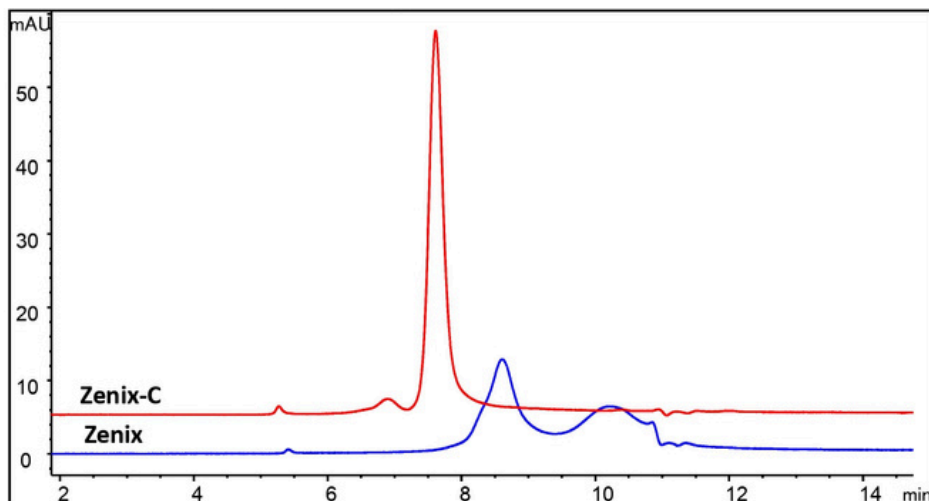
Detector: UV 280 nm;

Column temperature: 25 $^{\circ}$ C

Injection volume: 10 μ L;

Samples: Herceptin-lysine drug conjugate 2.05 mg/m

Antibody drug conjugate Analysis on Zenix-C 300

Column: Zenix SEC-300 (3 μ m, 300 \AA , 7.8 x 300 mm)

Mobile phase: 150 mM phosphate buffer, pH 7.0;

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
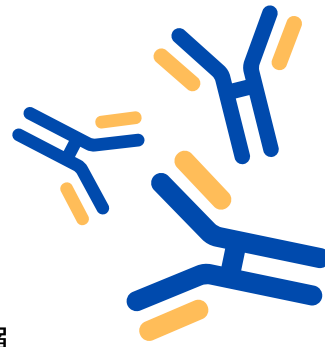
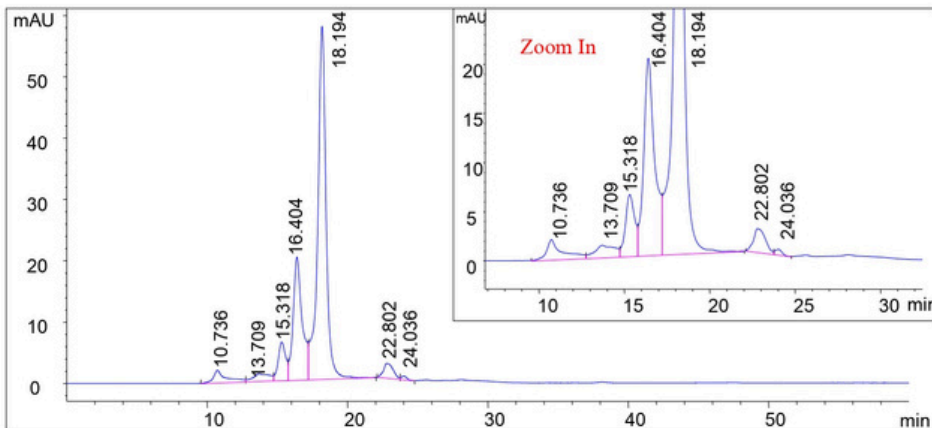
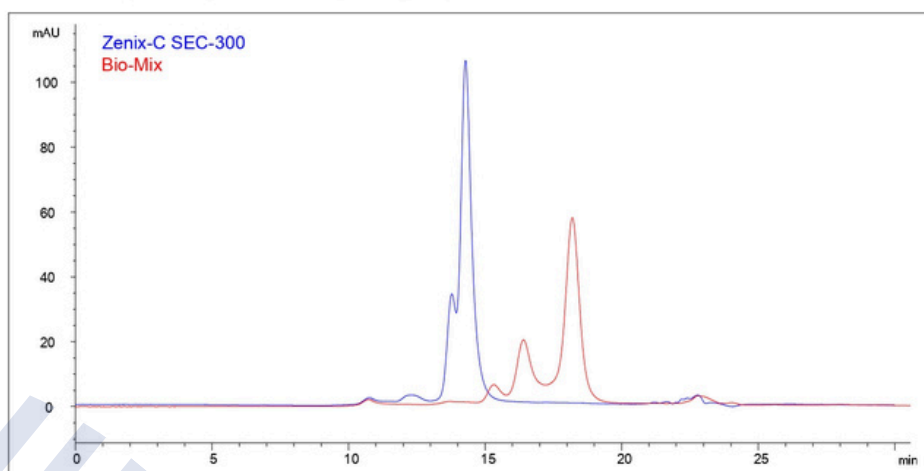
Operating Parameters:	
Operating temperature:	4-30°C
pH stability:	1-14
Chemical stability:	Resistant to aqueous solutions and most solvents used in liquid chromatography. Not resistant to acetone, ketones, chlorinated hydrocarbons, aliphatic esters, phenol, >10% NaOH, > 10% HCl, > 5% acetic acid, or strong mineral acid.
Materials:	
Glass column	Borosilicate glass
Endpiece	PTFE
Frit (bed support)	PE
O-Ring	FKM/FPM
Adjusting nut and retaining cap	Acetal
Connection cap and fitting nuts	Glass-filled polypropylene
Operating pressure:	
ID-6.6mm	900 psi (60 bar)
ID-10mm	600 psi (40 bar)
ID-15mm	300 psi (20 bar)
ID-25mm	150 psi (10 bar)
ID-35mm	150 psi (10 bar)
ID-50mm	100 psi (6.7 bar)

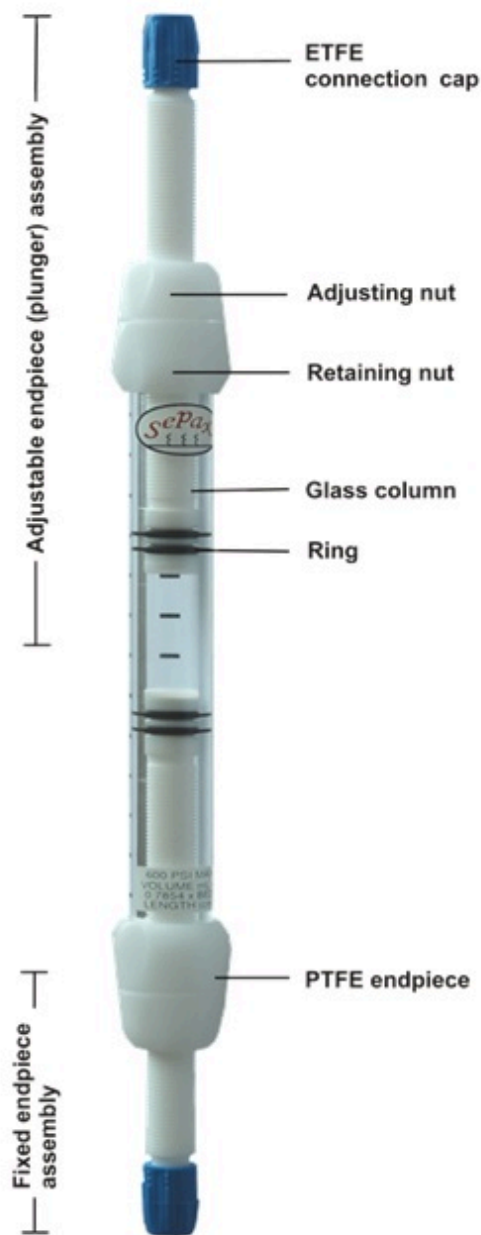3 configurations to choose from

AF: One fixed endpiece and one adjustable endpiece

AA: Two adjustable endpieces

FF: Non-adjustable with two fixed endpieces