# YOUR SEPARATION PARTNER

*Sepax Technologies*

Biopharma expert.

LINE Official Account

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@analab

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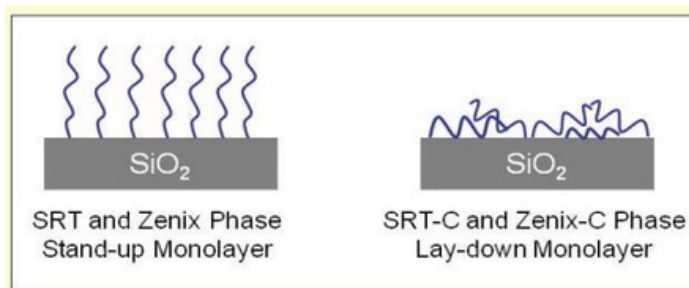
Ex: **hydrophobic proteins** like insulin, membrane protein, monoclonal antibodies derivatized with polymer branches, e.g. polypeptide, PEG.

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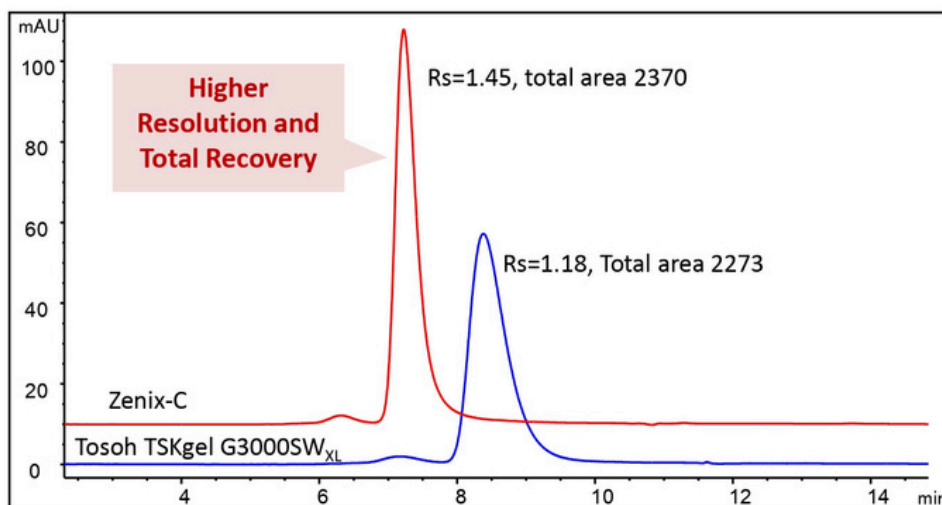
### Herceptin-Lysine Drug Conjugate Analysis

#### Zenix-C SEC-300 vs. Tosoh TSKgel G3000SWXL

**Column:** Zenix SEC-300 ( 3  $\mu$ m, 300  $\text{\AA}$ , 7.8 x 300 mm)

**Mobile phase:** 150 mM phosphate buffer, pH 7.0;

**Flow rate:** 1 mL/min;


**Detector:** UV 280 nm;

**Column temperature:** 25  $^{\circ}$ C

**Injection volume:** 10  $\mu$ L;

**Samples:** Herceptin-lysine drug conjugate 2.05 mg/m

### Antibody drug conjugate Analysis on Zenix-C 300

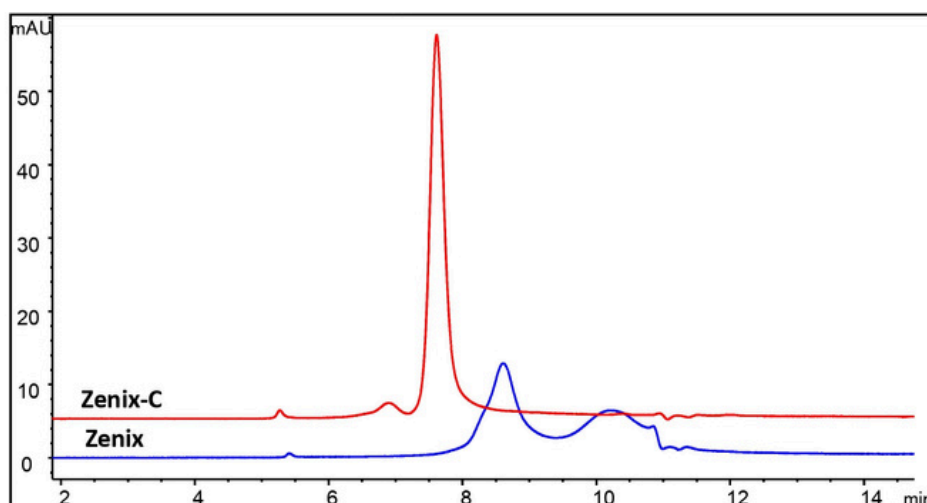**Column:** Zenix SEC-300 ( 3  $\mu$ m, 300  $\text{\AA}$ , 7.8 x 300 mm)

**Mobile phase:** 150 mM phosphate buffer, pH 7.0;

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
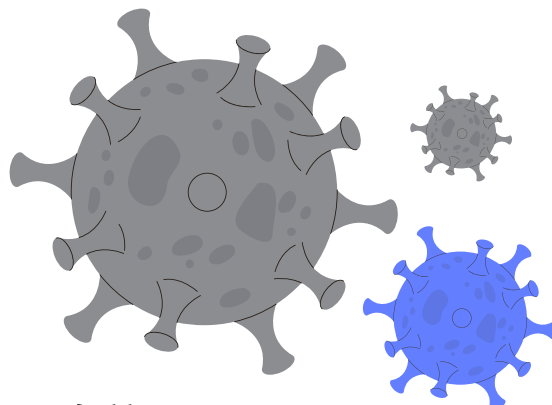
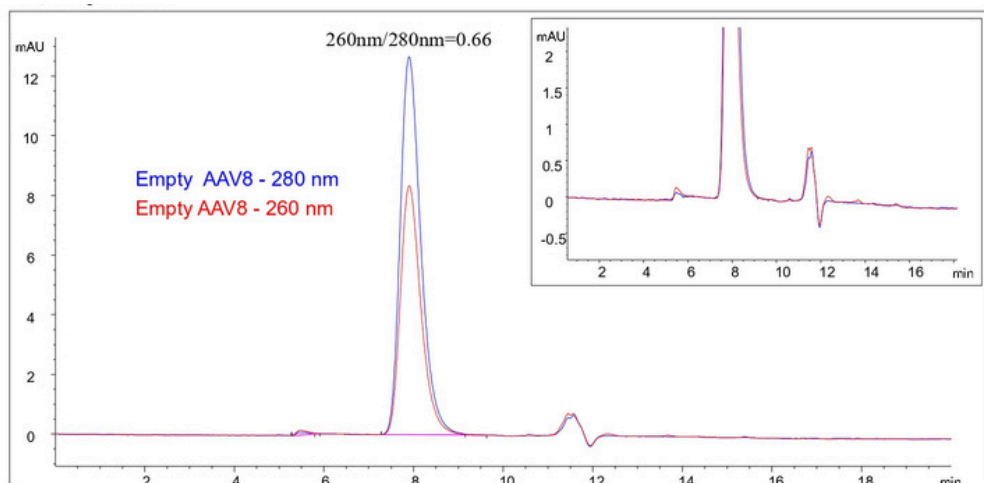
### AAV8 Pore Size Comparison

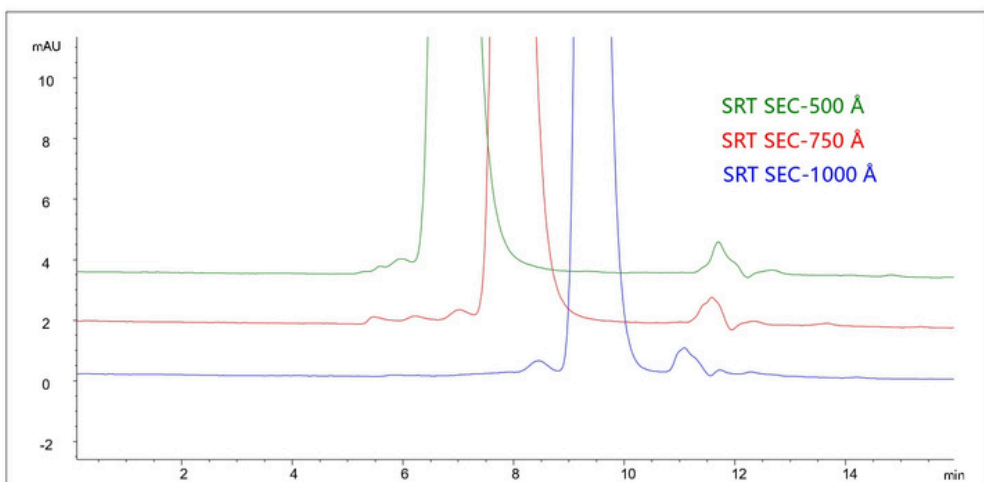**Column:** SRT SEC-750 Å, 5 µm, 4.6 x 300 mm;  
**Mobile Phase:** 1.8X DPBS with 10% Ethanol;  
**Flow Rate:** 0.35mL/min;  
**Column Temperature:** 25 °C  
**Injection Volume:** 10 µL;  
**Detector:** UV 260nm, 280nm;  
**Sample:** AAV8

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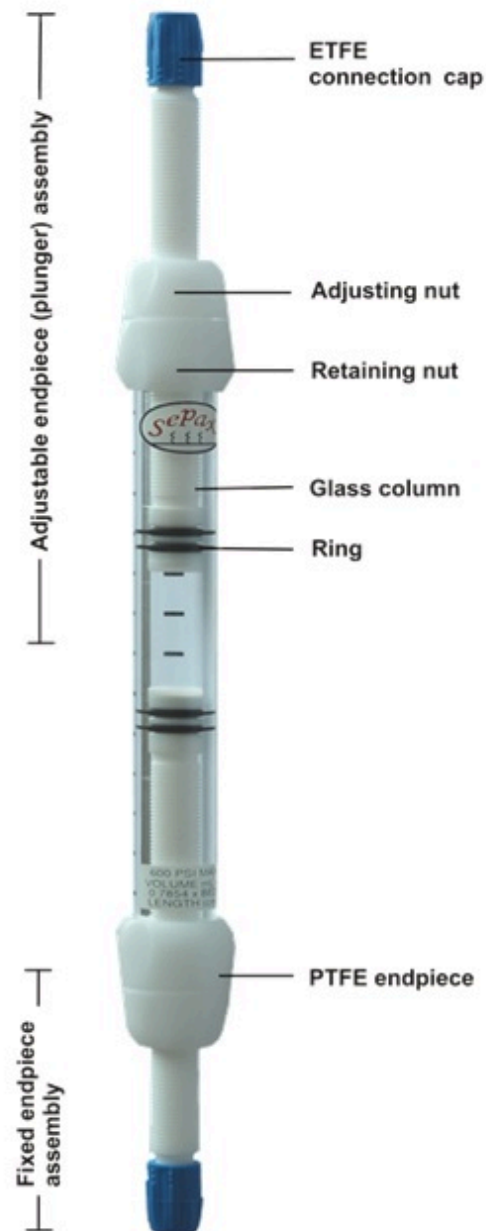
Operating Parameters:	
Operating temperature:	4-30°C
pH stability:	1-14
Chemical stability:	Resistant to aqueous solutions and most solvents used in liquid chromatography. Not resistant to acetone, ketones, chlorinated hydrocarbons, aliphatic esters, phenol, >10% NaOH, > 10% HCl, > 5% acetic acid, or strong mineral acid.
Materials:	
Glass column	Borosilicate glass
Endpiece	PTFE
Frit (bed support)	PE
O-Ring	FKM/FPM
Adjusting nut and retaining cap	Acetal
Connection cap and fitting nuts	Glass-filled polypropylene
Operating pressure:	
ID-6.6mm	900 psi (60 bar)
ID-10mm	600 psi (40 bar)
ID-15mm	300 psi (20 bar)
ID-25mm	150 psi (10 bar)
ID-35mm	150 psi (10 bar)
ID-50mm	100 psi (6.7 bar)


### 3 configurations to choose from

AF: One fixed endpiece and one adjustable endpiece

AA: Two adjustable endpieces

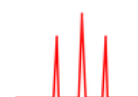FF: Non-adjustable with two fixed endpieces

立行科技有限公司  
Analab Corporation